

QUANTOFIX[®] Total iron 100

en

Pack content:

1 aluminium container with 100 test strips

Measuring range:

2 – 100 mg/L Fe²⁺ / Fe³⁺

Color gradation:

0 · 2 · 5 · 10 · 25 · 50 · 100 mg/L Fe²⁺ / Fe³⁺

General indications:

Remove only as many test strips as required. Close the container firmly immediately after removing the strips. Do not touch the test field.

Hazard warnings:

This test does not contain hazardous substances that must be labelled.

Instructions for use:

1. Dip the test strip into the test solution (pH 1 – 7) for 1 second.
2. Shake off excess liquid.
3. Wait 60 seconds.
4. Compare the test field with the color scale. If iron ions are present, the test field turns blue.

Interferences:

The coloration produced by some heavy metals in higher concentrations can be eliminated by adding cadmium powder.

Additional information:

The test strip container stopper contains a non-toxic drying agent. If swallowed, drink plenty of water.





Disposal:







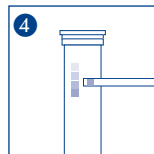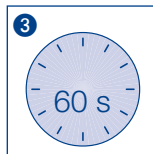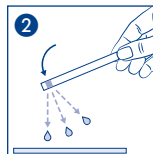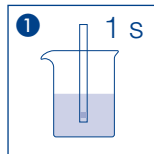


Information regarding disposal can be found in safety data sheet. You can download the SDS from www.mn-net.com/SDS.

Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4 – 30 °C).

If correctly stored, the test strips may be used until the use-by-date printed on the packaging.



 Artikelnummer / Item number / Référence produit / Referencia / Artikelnummer / Numero di catalogo / Numer artykułu	 Chargencode / Batch identification / Número de lot / Código de lote / Produktenummer / Codice del lotto / Numer partii
 Achtung! / Warning! / Attention! / ¡Atención! / Let op! / Attenzione! / Uwaga!	 Verwendbar bis / Use by / À utiliser avant / Fecha de caducidad / Te gebruiken tot / Utilizzare entro (anno / mese) / Przydatność do użycia
 Temperaturbegrenzung / Permitted storage temperature range / Limites de température / Limites de temperatura / Temperatuurbegrenzing / Limite de tem- peratura / Ograniczenie temperatury	 Trocken aufbewahren / Store in a dry place / Conserver au sec / Mantener seco / Op een droge plaats bewaren / Manteneren asciutto / Przechowywać w suchym miejscu
 Gebrauchsanweisung beachten / Please read instructions for use / Respecter les instructions d'utilisation / Obsérvese las instrucciones de uso / Lees de bijsluiters / Consultare le istruzioni per l'uso / Przestrzegać instrukcję użycia	 Packung geschlossen halten / Keep container closed / Refermer la boîte / Mantenga el envase cerrado / Verpak- king gesloten houden / Conservare la confezione chiusa / Przechowywać pojemnik szczelnie zamknięty
 Ausreichend für <n> Prüfungen / Contains sufficient for <n> tests / Contenu suffisant pour <n> tests / Contenido suficiente para <n> tests / Voldoende voor <n> tests / Contenuto sufficiente per <n> test / Wystarczająco dla kontroli <n>	